




Duhlwiesen 32, 55413 Weiler bei Bingen am Rhein - Germany
Tel. +49-6721-9886710, Fax. +49-6721-9886719

Technical Note

TN0002

TDAT Problem Solutions Guide

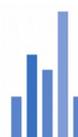

Titel

TDAT Problem Solutions Guide

Editor

Oliver Mueller / Parts Engineer

Version


1.00

Updated

19.11.2009

Instruments

MassLynx Systems

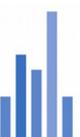

Contents

1 TDAT Introduction.....4

2 TDAT Acquisition Cards.....5

3 How to diagnose a TDAT card.....7

4 Introduction to EngCon.....8

5 Additional information.....11

6 How to build a TDAT Cable.....12

7 Error Codes.....13



1 TDAT Introduction

This guide is dedicated to all users, who still have a Micromass TDAT instrument in use. Actually there is not much knowledge left from the good old Micromass times and the old engineers might be able to remember it. This guide is a collection of information what Waters provided for their customers. Have a look in the manuals that come with a new TDAT PC or what you find in the Masslynx 3.5 knowledge base help. All pictures are made by ourselves and all information are given with no claim to be complete. For your reference, have a look in the Waters document 715001758 Rev. A.

Since the early Trio-family made by VG-Organic, TDAT was a standard interface and used for instrument control and data acquisition. The TDAT card in the Masslynx PC communicates with the MPU or TPC board in the Mass Spectrometer. It uses a twisted pair serial communication line. The name TDAT is derived from Trio-Data-Acquisition. Later it was used by Micromass and (Finnigan). Since the year 2000 the TDAT control was substituted by the embedded PC acquisition system (EPCAS).

There are several disadvantages with TDAT, for example limited cable length between the instrument and the control PC (max. 5 meters). The placement of the cable could cause communication problems especially when it is close to strong magnetic fields like roughing pump motors or transformers. The other limitation with TDAT based systems is the availability of the microprocessors. Most of them were produced in the late 80's. Each instrument required two controllers. One is the SPU or Scan Processing Unit. It is responsible for instrument communications and control. The other is called the PPU or Peak Processing Unit that is responsible for data acquisition and the tune page display. The two processors were later combined on one board called the MPU or Multiple Processor Unit. This was only used on the Platform I and II. It was then updated to the TPC or Transputer PC board that is used in all newer products.


Some TDAT instruments are supported in Masslynx 4.1 but only with use of a PCI TDAT card.

For more information please ask your local Waters representative. For some Quattro LCs and Ultimas upgrade kits are available.

2 TDAT Acquisition Cards

ISA TDAT Card

TN0002 TDAT	
-------------	--

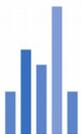




This picture shows the TDAT ISA card. Notice the two sliders for the IRQ (interrupt request) and DMA (direct memory access) settings. By default set IRQ 5, DMA 1 and the Jumper to 2.5 MHz. You can use the windows device manager to configure the card's resources. A working card will not be listed in the installed hardware tab. Check in the Device Manager, if a IRQ and DMA number was assigned. If you need to change the settings you need to change the parameters in the registry, run regedit, go to Hkeylocal machine/system/currentcontrolset/services and search for the TDAT entry and alter the parameter tab. After rebooting the PC, windows will try to start the tdat.sys driver with the new settings. If your card will fail again, even though in a working computer, you might need to replace the large foil capacitor.

PCI TDAT Card



The PCI TDAT card was made for use in modern PCs. The bad thing is, that this card is not a true plug and play card and you need special software change notes. First you install Windows with standard PC drivers. You can change the settings in the device manager, click on the computer icon and change the driver from ACPI PC to

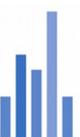


Standard PC. Windows needs to be rebooted and will reinstall all drivers. Then you should install Masslynx and the required software change note (SCN). Shut down the PC and insert the PCI TDAT card. After the new start Windows finds a new device.

ML4.0 first release needs SCN 433

ML4.0 SP2 needs SCN 485

ML4.0 SP4 needs SCN 508

The SCN includes the driver that is copied to the Masslynx folder. With the driver comes a small program tdattest.exe that is for TDAT card diagnostics.

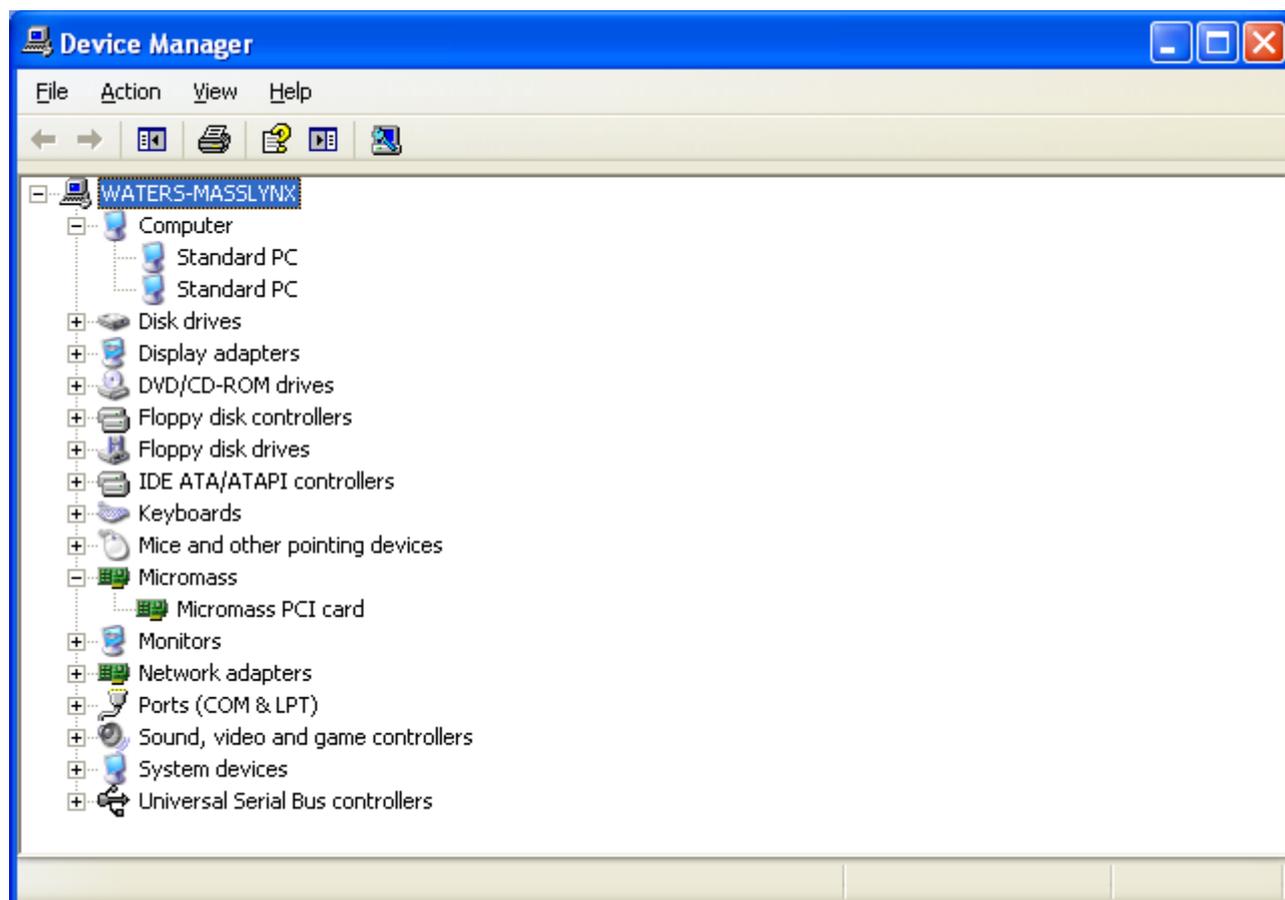

Tip: If you need to buy a PCI TDAT card, Waters charges thousands of euros for it. The card itself is no longer available as a single card, the part number disappeared and you need to order the upgrade kit. This Kit is 3 times more expensive as the card alone was, but it contains a CD with the SCN and a short manual. The best idea is to buy a complete TDAT PC from Waters for almost the same price as the kit.

3 How to diagnose a TDAT card

You need a TDAT loopback connector that plugs into the 15 pin port of your TDAT card.





The pictures show this connector, it is made very simple.

Pinout for 15 Way D-type loopback connector

Link Pin 1 to Pin 10

Link Pin 2 to Pin 9

Link Pin 3 to Pin 8

Link Pin 4 to Pin 7

Link Pin 5 to Pin 6

With the connector installed, start TdatTest.exe

```
C:\Masslynx\TdatTest.exe
Failed to initialise Transputer (status =0)
Failed to identify transputer (got cstatus = 81,ctype = 0)?

-----
PCI/ISA IDAT Tests - Main Menu
L ..Loopback using program I/O *needs loopback connector
D ..Loopback using DMA *needs loopback connector
T ..Loop on initialise & identify Transputer
R ..Read/Write Transputer memory
E ..Exit to Windows


Loopback test ...
Loopback test done...Pass 1144
```

Here you can choose L or D to test the card. It will increment the Passes until you hit a key. To diagnose a card it might be necessary to run the test a couple of hours.

4 Introduction to EngCon




EngCon (engineer control) is a small program located in the Masslynx folder to control the mass spectrometer directly. It is only available on TDAT systems.

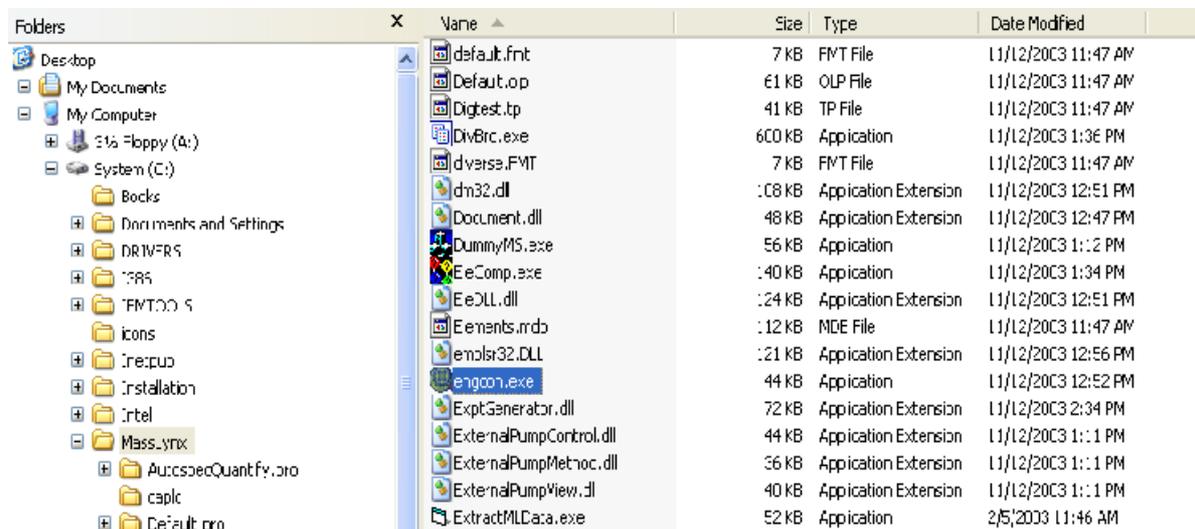


Folders	Name	Size	Type	Date Modified
Desktop	default.fmt	7 KB	FMT File	11/12/2003 11:47 AM
My Documents	Default.op	61 KB	OLP File	11/12/2003 11:47 AM
My Computer	Digitest.tp	41 KB	TP File	11/12/2003 11:47 AM
3 1/2 Floppy (A:)	DivBrc.exe	600 KB	Application	11/12/2003 1:36 PM
System (C:)	dverse.FMT	7 KB	FMT File	11/12/2003 11:47 AM
Books	dm32.dll	108 KB	Application Extension	11/12/2003 12:51 PM
Documents and Settings	Document.dll	48 KB	Application Extension	11/12/2003 12:47 PM
DRIVERS	DummyMS.exe	56 KB	Application	11/12/2003 1:12 PM
RAM	EeComp.exe	140 KB	Application	11/12/2003 1:34 PM
FMT000	EeDLL.dll	124 KB	Application Extension	11/12/2003 12:51 PM
icons	Elements.mdb	12 KB	MDE File	11/12/2003 11:47 AM
irecpu0	emolr32.DLL	121 KB	Application Extension	11/12/2003 12:56 PM
Installation	engcon.exe	44 KB	Application	11/12/2003 12:52 PM
Intel	ExptGenerator.dll	72 KB	Application Extension	11/12/2003 2:34 PM
Masslynx	ExternalPumpControl.dll	44 KB	Application Extension	11/12/2003 1:11 PM
AucSpecQuantfy.pro	ExternalPumpMetroc.dll	56 KB	Application Extension	11/12/2003 1:11 PM
caplc	ExternalPumpView.dll	40 KB	Application Extension	11/12/2003 1:11 PM
Default.pro	ExtractMLData.exe	52 KB	Application	2/5/2003 11:46 AM

If you want to start EngCon, you should close Masslynx because only one program at a time can have access to the TDAT control. Double click on engcon.exe opens this program and you can open a transputer file for your instrument. If you have an Ultima, choose Ultimam.tp. For Platform or ZMD choose a file named platicz.tp. ZMDs show an instrum.cfg error (02) message. That is because the *.tp file was developed for the LCZ. Anyway, it will work.

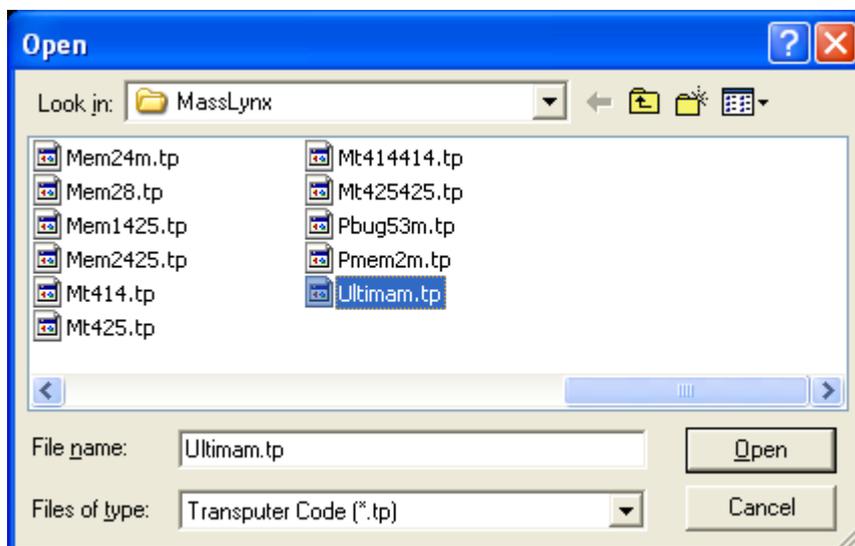

After clicking “open” you will receive a window like the one below. EngCon does a couple of self tests and when it is ready, hit the Escape (Esc) key.

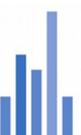

The image shows two screenshots of the 'Ultimam.tp - NT EngCon' software interface. The top screenshot displays the 'SPU : Setup - STARTED' process, listing various initialization steps such as 'init.a4d.board', 'init.csrs.etc', and 'activating PPU', all of which are marked as 'FINISHED'. The bottom screenshot shows the 'MISC Page GAS' configuration screen, which includes a copyright notice for 'Masslynx QUANTRO ULTIMA v3.5_1' and a table of instrument settings.

```

SPU : Setup - STARTED
TracePackage Init: TRACE.ENABLED=FALSE, TraceLevelCurrent=TraceLevelOverride
SPU : Setup - init.a4d.board - FINISHED
SPU : Setup - init.csrs.etc - FINISHED
SPU : Setup - getting imr ..... imr reported as 2000
SPU : Setup - get.config.info - FINISHED
SPU : Setup - activating PPU - FINISHED
SPU : Setup - performing power.up.self.test
No GC detected - FINISHED
SPU : Setup - Initialising global variables - FINISHED
SPU : Setup - IEnergy Ramping Disabled ... Initialising DACs - FINISHED
SPU : Setup - FINISHED

```

```


MISC Page GAS
Masslynx QUANTRO ULTIMA v3.5_1
Copyright 1999 Micromass Ltd.
All Rights Reserved
4190 AMU QUAD

[OP] OPERATE OFF [TO] TURBOS ON OFF
[CC] COLLISION CELL GAS OFF [PC] PUMP COLLISION CELL OFF
[E2] APCI/ESI GAS OFF [AR] AutoRanging a4d ON
[IL] INJECT ON [PR] PRESSURE REQUESTED OFF

<ESC> to select another page or <CTRL> X to exit
SRC / ION / SCAN / ACQU / MISC / MTHD / RDB / INLT / ASC / MS2
>on

```

If your MS is in vent state, you can navigate by using the arrow keys. Go to misc/gas and move to the highlighted [TO] Turbos ON / OFF. You will notice the prompt > in

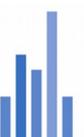

the command line. To start the roughing pump and the turbo pumps type on and hit enter, type off to vent the MS. Under RDB you find the readback pages for the turbo speed. Use the arrow keys for navigation or type rd1 for page1, rd2 for page 2, and so on.

What is the use for this program?

You can use it to check voltages, align the RF-Generator, do other instrument tests. There is the possibility to do MS-Scans. This is only used for board setup procedures, not for data acquisitions like you know from the tune page.

Most used commands

G [RETURN]	to start scanning
H [RETURN]	to stop scanning
I [RETURN]	to go to tests
T [RETURN]	to start Mass Program
V [RETURN]	shows instrument related variables

5 Additional information


During the Masslynx installation you will be asked what mass range your spectrometer has, or if you have the cone gas option installed and so on. All this settings are stored in the Instrum.cfg file. The most important one is the ZZ variable that determines the mass range. For a Qultima the file is like the one below.

```
MD=16
ZZ=4190
RO=0
RL=400
RR=250
ZS=1
R1=0
R2=0
```

The type of your mass spectrometer is stored in the AI.ini file.

```
[Boot]
BootFile=ULTIMAM
```

```
[Options]
PauseBoot=0
IconStart=1
```

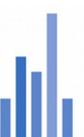

Please make sure that the files match your instrument.

6 How to build a TDAT Cable

First you need to buy a high quality shielded cable to keep the signal noise low. The maximum cable length should not exceed 5 meters.

Note: The cable pinout is numbered from the view of the connector solder site.

MS-Connector		TDAT-Card-Connector	
Link Pin	1	to Pin	8
Link Pin	2	to Pin	7
Link Pin	3	to Pin	6
Link Pin	4	to Pin	5
Link Pin	7	to Pin	2
Link Pin	8	to Pin	1
Link Pin	9	to Pin	15
Link Pin	10	to Pin	14
Link Pin	12+13	to Pin	11+12+cable shield
Link Pin	15	to Pin	9



7 Error Codes

The transputer card in the mass spectrometer has two red multisegment displays. If everything is o.k. they show 0,0. In case of an error the instrument will display an error number.

Error number list

- 01 - Missing CAL/CFG file: Cannot find INSTRUM.CFG or a .CAL file
- 02 - Bad CAL/CFG file format: Bad INSTRUM.CFG or .CAL file format
- 03 - Bad GC/Autosampler reply: Bad or no reply from GC or Autosampler
- 04 - DuartIntBadCh: DUART interrupt received but wrong character
- 05 - DuartNoInt: DUART didn't generate interrupt (faulty chip?)
- 06 - DuartIntNoCh: DUART interrupt received but no data
- 07 - WrongEventDuart: DUART generated interrupt but got wrong event (
- 08 - EventFailDuart: DUART interrupt not detected (faulty event reg?)
- 09 - DuartLoopFailB: DUART local loopback failed port B (faulty chip)
- 0A - DuartLoopFailA: DUART local loopback failed port A (faulty chip)
- 0B - DuartNotThere: DUART not responding
- 0C - DuartSpuriousInt: Unsolicited interrupt from DUART
- 0D - TxServerFail: Cannot transmit to PC server
- 0E - RxServerFail: Timeout on waiting for PC data
- 0F - ServerResetFail: Server reset didn't work - server corrupted or in unknown state
- 10 - BadResetChecksum: PC data corrupted or bad communications



- 11 - Missing PPU.MPO file: Cannot find TMS chip code file
- 12 - Bad PPU.MPO file format: TMS code chip file has bad format
- 13 - CAL coefficient(s) out of range: Coefficient(s) out of range in a .CAL file
- 14 - Wrong code for scan board: Quattro/Bio-Q has been booted with wrong code for scan board fitted
- 15 - No scan board fitted: No scan board could be detected on a Quattro/Bio-Q
- 16 - Scan board detected failed: Quattro/Bio-Q scan board detection failed because of hardware fault
- 2A - Inter-scan overrun: Transputer too busy during inter-scan delay to start next scan on time
- 2B - Scan not finished: Hardware taking too long to do scan (hardware fault?)
- 2C - Unrecognised event: Unrecognised type of event interrupt recieved
- 2D - Can't open datafile: Could not open a file via server
- 2E - Error writing data: Error while writing data to file
- 2F - Data buffers full: Scanning too fast or cannot get enough access to disk
- 30 - Bad checksum:Corrupt server communications
- 31 - Bad record length: Corrupt server communications
- 32 - Bad server command:Corrupt server communications
- 33 - ASStatus message: Receeived status message from Autosampler
- 34 - Undefined Error Code
- 35 - ADC still busy: ADC conversion not completed in time allowed
- 36 - Timeout server comms: server and/or link failure
- 37 - PC disk full: PC disk full - acquisition halted
- 38 - Filelist disk error: Disk error during filelist processing
- 39 - Filelist format error: Filelist control file has bad format
- 3A - Filelist command error: Error in filelist command section
- 3B - Undefined Error Code
- 3C - Total communications failure: Server or link dead
- 3D and 3E Failed to identify sampling rate of the AutoRanged Analog channels (is the jumper fitted correctly)ed to EMC
- 3E - AutoA4DFifoFull AutoRanged FIFO has filled up due to irregular servicing (see table below for maximum scan times)

The following is based on me interpreting the code on the Quattro LC.



On the Quattro LC, 3F is a 'Calibration timeout error'. This occurs at the start of a scan, when the DACs (MS1, Ramps. MS2) are calibrated. The software clears some bits what it calls 'digital IO' and then waits 250 ms for the board to register that it has done what it was supposed to. If it fails to, then it registers the error.

The code mentions the AD760 data sheet, if that helps.

FF is a bit more tricky to suss out, but I've seen it a lot before now on perfectly happy systems. It appears to be a default status that turns up, particularly when you connect MassLynx or start acquisitions.

5* errors are related to all acquisitions and may be due to acquiring saturated beams, or too much data too quickly.

- 51 - Bad data flag from TMS: TMS program misbehaving
- 52 - FIFO 1 overrun: Acquiring too fast / TMS hung / Transputer slow
- 53 - No response from TMS: TMS program crashed or hardware fault
- 54 - Bad start flag from TMS: TMS program misbehaving
- 55 - Undefined Error Code
- 56 - FIFO 2 overrun: Acquiring too fast or too much data

6* errors are related to centroid scans and may be due to acquiring saturated beams, or too much data too quickly.

- 64 - Overflow in moment sums
- 65 - Overflow in average sums
- 66 - Overflow in filter calculation
- 67 - Overflow in derivative calculation
- 68 - Overflow in time calculation
- 69 - Untracked overflow
- 6A - Bad transputer scan code
- 6B - Accumulator overflow
- 6C - FIFO 1 overrun
- 6D - Bad transputer scan parameters



E* messages are related to the boot emproms fitted to Quattro I, Quattro II and Trio-200 instruments.

The following codes appear in sequence as the machine boots up and appear for less than five seconds, after which time MassLynx behaves normally.

E0 - EpromBootStart

The machine is starting to initialise using internal EPROMs to boot after a power on

E1 - EpromBootCSR

Initialising Vent/Standby/user IO ports

E2 - EpromBootSRC

Initialising SRC page as for ESP source

E3 - EpromBootSCN

Initialising SCN page to low masses

E4 - EpromBootMSC

Initialising miscellaneous dacs such as RF lens/Collision

Energy/Gas Cell to zero ramps

EB - EpromBootEnd

Finished initialisation

